

BetterBuffer

Sequence enhance/dilution buffer

Specially formulated to be used with ABI BigDye™ Mix (versions 2.0 and 3.0)

Protocol:

Mix **5** µl of **BetterBuffer** with **1** µl of ABI BigDye™ mix*

Add DNA and primers to a final volume of 15 µl

Sequence as normal

Spin out the excess dye with ethanol or 80% isopropanol

BetterBase

Sequence enhance/dilution buffer

Specially formulated to be used with ABI BigDye™ Mix (versions 1.1 and 3.1)

*other ratios may be preferred

BigDye is a trademark of PE Applied Biosystems

Store at -20°C

For Research Only